

Experiment 1: Introduction to Labview 8.0

(tbc 1/7/2007, 1/13/2009,1/13/2011)

Objective: To learn how to build virtual instrument panels in Labview 8.0

Tasks:

1. Build the Labview VI shown in the tutorial given in Appendix A.
2. Modify the VI such that:
 - a) The formula is now given by
$$x = A \sin(\omega_1 t) + B \cos(\omega_2 t)$$
 - b) The input nodes are now: amplitudes A and B , frequencies ω_1 and ω_2 , and elapsed time t . This means you need to include three more controls that allow the user to adjust the first four variables.
 - c) Instead of sliders, use **[Numeric Ctrl]** controls. This means you do not need to attach a numeric indicators to the “Numeric Ctrl” controls. Further, these allow the user to input values directly. Change the properties of the “Numeric Ctrl” controls to match Table A. (see Figure A.)

Table A.

Numeric Ctrls: A and B	Minimum	0
	Maximum	0.5
	Increment	0.01
Numeric Ctrls: w1and w2	Minimum	0
	Maximum	10
	Increment	0.01

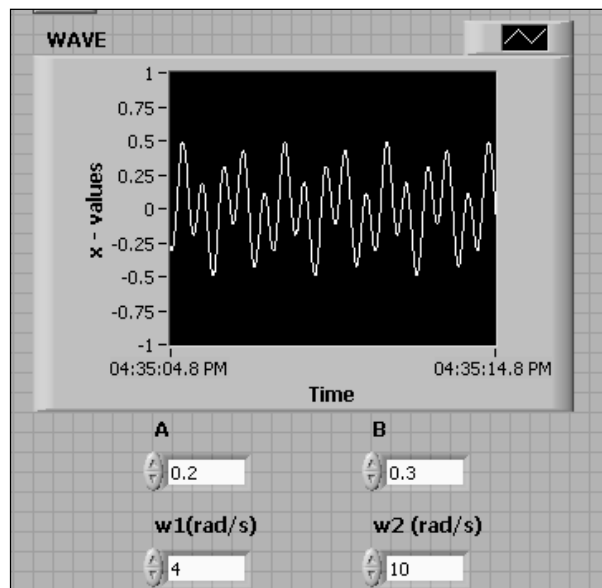


Figure A. Modified Front Panel

Appendix A: Tutorial for Labview 8.0

I. Initializing a blank virtual instrument.

1. Log on to **Labview 8.0** and open a **Blank VI**. (see Figure 1.)

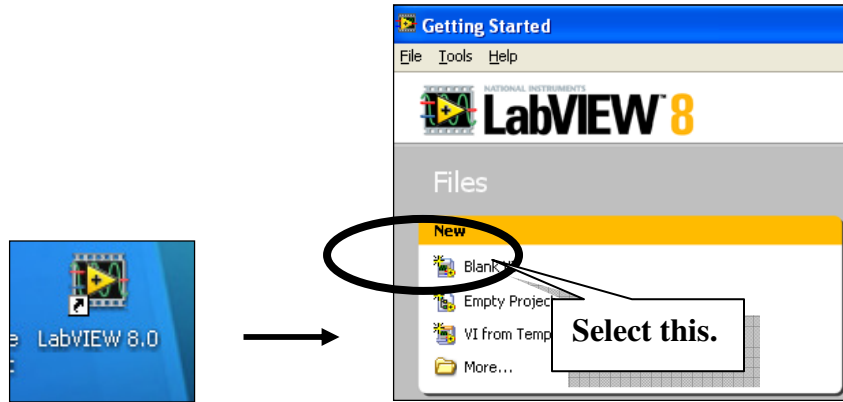


Figure 1. Opening a blank virtual instrument (VI).

2. Two windows will appear: a) the **Front Panel** and b) the **Block Diagram**. (You could tile both windows side-by-side by selecting [**Window**]→[**Tile Left and Right**] menu item as shown in Figure 2).

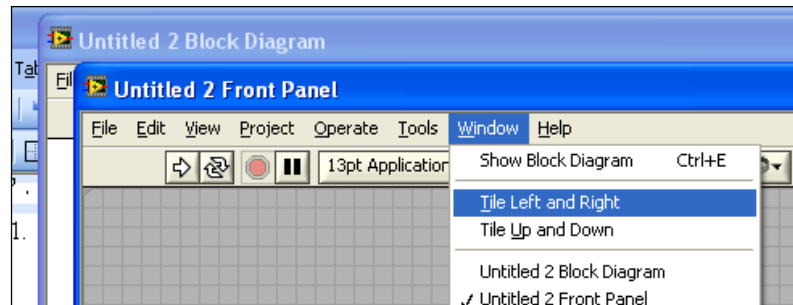


Figure 2. The “Front Panel” and “Block Diagram” windows.

Remarks:

- i) The “Front Panel” will contain the user interface such as knobs, buttons, indicators and graphs while the “Block Diagram” window will contain the inner-circuitry (including objects, wirings and other virtual instruments), logic structures (including programming blocks) and timing commands.
- ii) If either the “Front Panel” or the “Block Diagram” window is blocked by another window (e.g. when either window is maximized), one can toggle the focus between the two windows by pressing [**CTRL**]-[**E**].

II. Building a VI. (Sliders, Functions and Charts)

Description: We will now build a virtual instrument (VI) project that produces a sinusoidal signal whose frequency will be adjusted by a slider. We will also monitor the signal using a strip chart and save the results in a text file.

Procedure:

1. Access the Functions-palette.

[Right-click] on the mouse cursor while inside the **block-diagram** window, then “tack”-on the **Functions**-palette (see Figure 3).

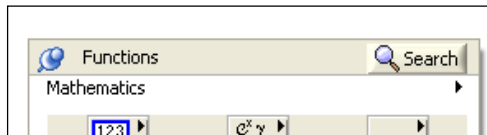


Figure 3. Tack-on by clicking on the the push-pin.

2. Create a While-Loop box.

By clicking on the dark triangles in the Functions palette, one can expand or reduce different icon-subdirectories of available functions.

Expand to the **[Express]→[Exec Control]** subdirectory, then click-on the **[While-loop]** icon and go to the block-diagram window to create an area for the while-loop function as shown in Figure 4.

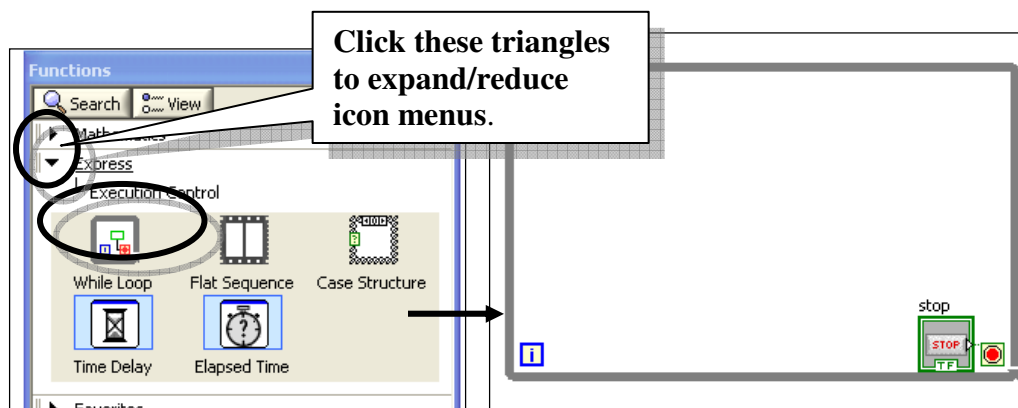


Figure 5. Click the while-loop icon, then drag an area in the “Block Diagram” window.

Remarks:

- i) A **[STOP]** button will also appear in the **front-panel** window.
- ii) You can have several Function palettes open at the same time.
- iii) You could select or de-select “Function Groups” (e.g. Mathematics, Express, etc.) by checking or un-checking the items under the **[View]→[Always Visible Categories]** submenu, as shown in Figure 6.

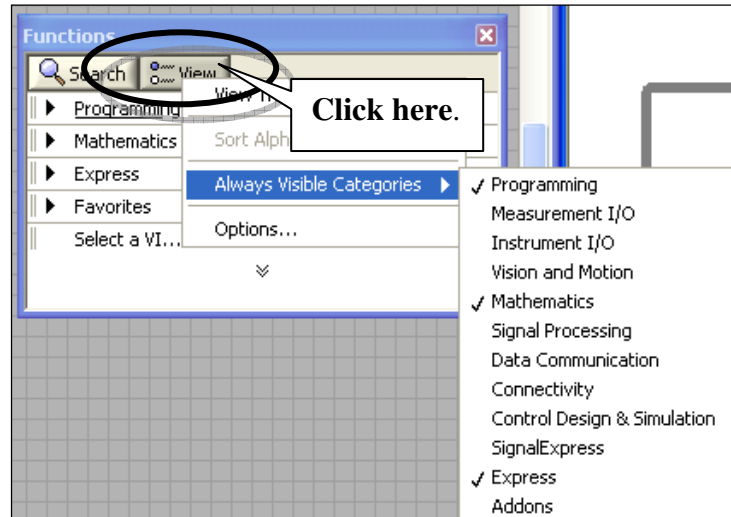


Figure 6. Selecting and deselecting function groups.

NOTE: For the items below, make sure that all the objects created are inside the **[while-loop]** box.

3. **Build a Formula Node.**

In the Functions palette, expand the **[Mathematics]→[Scripts & Formulas]** icon subdirectory, then click on the **[Formula Node]** icon. In the “Block-diagram” window, drag an area then enter the formula: **$x = \sin(w * t);$** as shown in Figure 7. (Remember to include the semicolon).

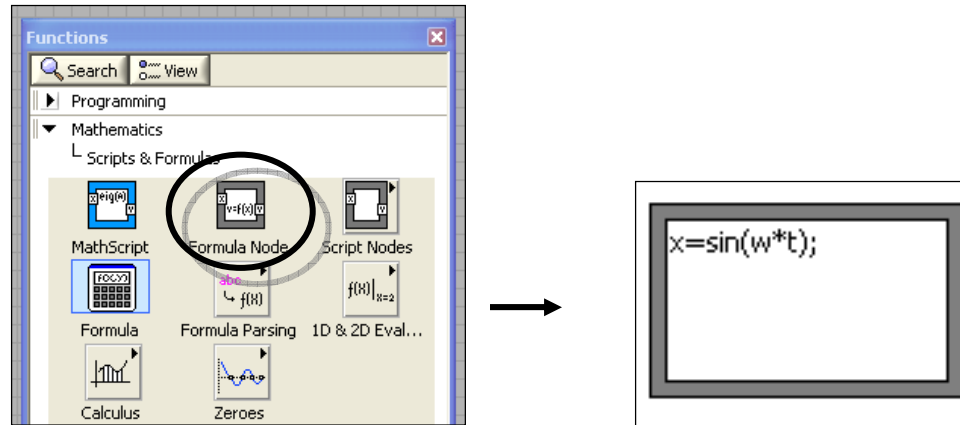


Figure 7. Introduce a Formula node block.

Move the cursor to hover around the left edge of the “Formula-node” block, then **[Right-Click]** and select **[Add Input]**. A small box should appear. Name it: “**w**”. Add another input node and name it: “**t**”. Similarly, hover around the right edge of the “Formula-node” block, **[Right-Click]** and select **[Add Output]**, then name it: “**x**”. (The output nodes will have a thicker boundary, see Figure 8).

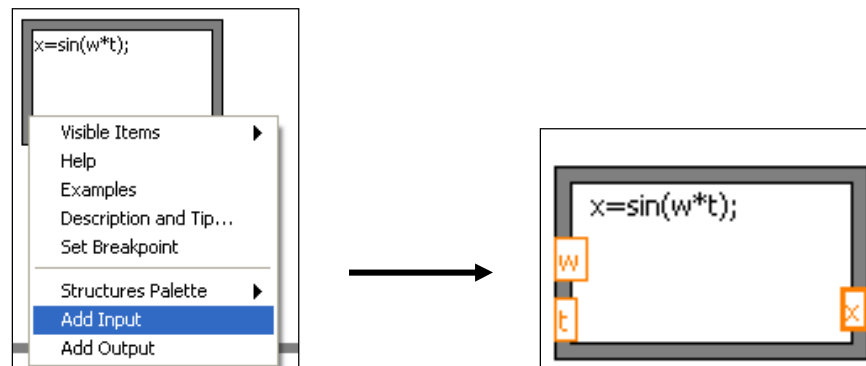


Figure 8. Adding input and output variables.

4. **Insert the “Elapsed Time” block then connect it with the “t” input.**

From the Functions palette, expand the **[Programming]→[Timing]** icon subdirectory and select the **[Elapsed Time]** icon. Drop the object inside the “Block-diagram” window. A “**Configure Elapsed Time...**” window should appear. Disable the **[Automatically reset...]** checkbox, then click the **[OK]** button. (See Figure 9.)

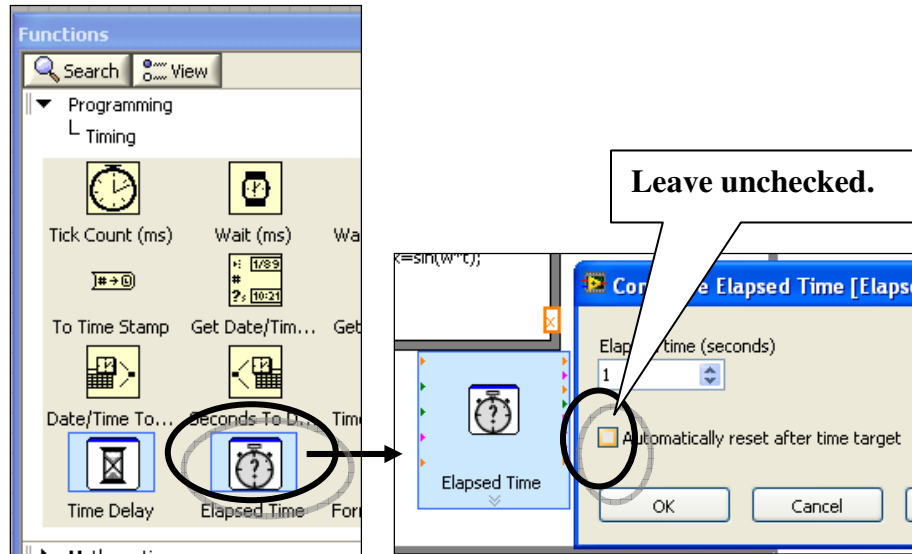


Figure 9. Insert an “Elapsed Time” block.

Let the cursor hover around the “**Elapsed Time (s)**” selection until the cursor changes to a “wire-spooling” cursor. Next, **[Click-drag]** the spool until it hovers near the “**t**” input node of the “Formula-node” block and then release. A connecting wire should then appear. You may need to resize the “While-loop” box and/or move the “Formula-node” block to make room for the desired wiring between the “Elapsed Time” block and the “Formula-node” block. (See Figure 10).

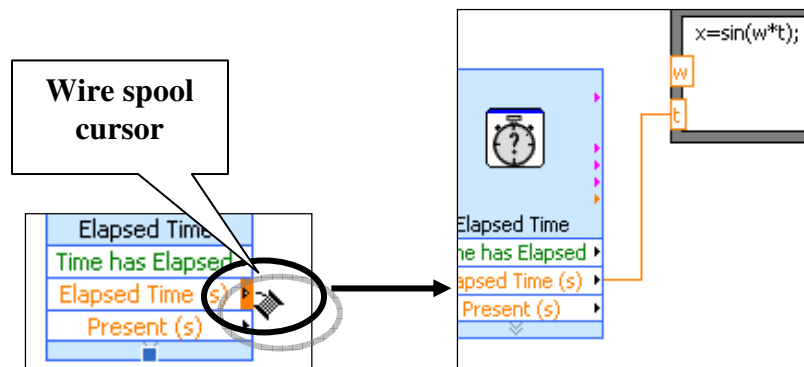


Figure 10. Wiring the “elapsed time” values to the “t” input node.

Remarks:

- i) To save space, you could **[Right-click]** the “Elapsed time” block and select the **[View as Icon]** menu item. This should reduce the block to be an icon (See Figure 11.)

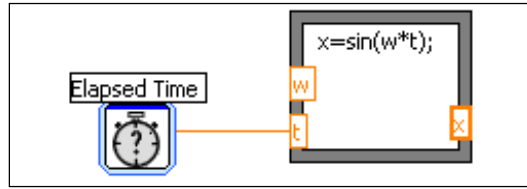


Figure 11. “Elapsed Time” block reduced into an icon.

- ii) When the cursor is in the “pointer” mode, a **[Click-drag]** could be used to move or remove blocks and wires.

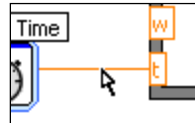


Figure 12. “Pointer” cursors.

- iii) Frequently, a wire may become “broken”. This is usually marked by an “X” on the broken wires as shown in Figure 13. This means that the broken wire that has one unattached end. To remove all broken-wires in the “Block-diagram” window, press **[CTRL-B]**.

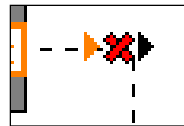


Figure 13. A broken wire.

5. Add a horizontal “Slider” control.

[Right-Click] at a region in the “Front Panel” window and then **[Click]** the push pin to access the “Controls palette”.

Expand in the “Controls palette”, the **[Express]→[Numeric Controls]** icon subdirectory, then select the horizontal **[Pointer Slide]** icon and drop it inside the “Front Panel” window. **[Double-Click]** the “Slider” label and rename it as as “Frequency (rad/s)”. (See Figure 14.)

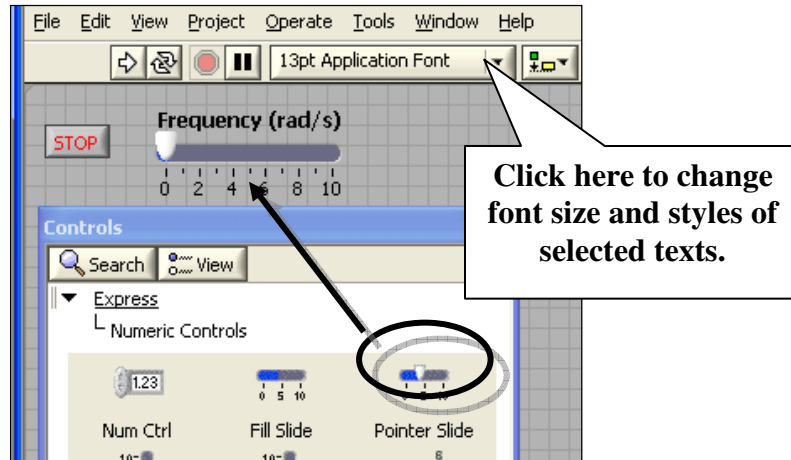


Figure 14. Setting a horizontal slider.

Next, go to the “Block Diagram” window and make sure that the “Frequency (Slider)” block is inside the “While-loop” box. (If not, drag it inside.)

Move the cursor around the small output triangle node at the right edge of the “Frequency (Slider)” block until the “wire-spool” cursor appears, then drag a wire to connect the “Frequency (Slider)” block to the “**w**” input node of the “Formula node” block. (See Figure 15.)

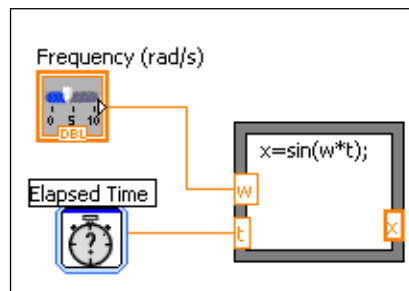


Figure 15. Connecting the output of “Slider” block to one of the inputs of “Formula node” block.

6. Add a “Chart” control.

[Click] at a region in the “Front Panel” window and access the “Controls palette”. Under the **[Express]→[Graph Indicators]** icon subdirectory, select the **[Chart]** icon then drop a “Chart” control in the “Front Panel” window. Rename the “Chart” control to “SINE WAVE”. (See Figure 16).

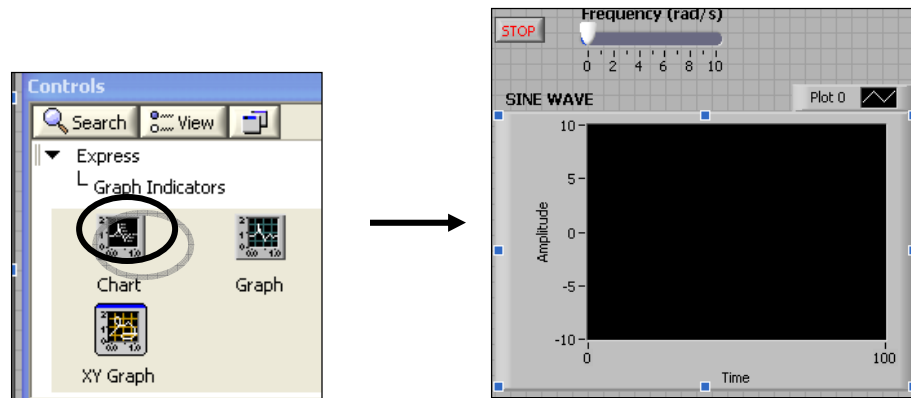


Figure 16. Adding a “Chart” control inside the “Front Panel” window.

Next, go to the “Block Diagram” window and make sure that the “SINE WAVE (Chart)” block is inside the “While-loop” box. (If not, drag it inside.)

Move the cursor around the small input triangle node at the left edge of the “SINE WAVE (Chart)” block until the “wire-spool” cursor appears, then drag a wire to connect the “SINE WAVE (Chart)” block to the “**x**” output node of the “Formula node” block. (See Figure 17)

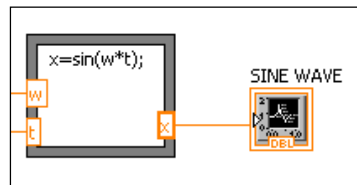


Figure 17. Connect the “**x**” output node to the “SINE WAVE” block.

7. Add a “Wait Until Next Multiple ms” block.

Access the “Function palette” and expand the **[Programming]→[Timing]** icon subdirectory. Select the **[Wait Until ...]** icon, then drag it into the “Block Diagram” window. (See Figure 18).

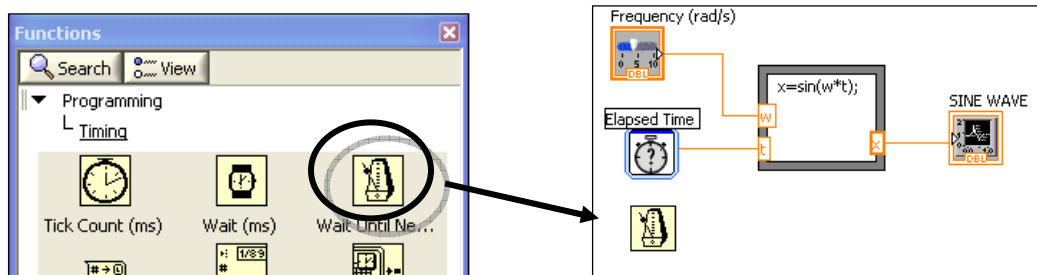


Figure 18. Add a “Wait ...” block into the “Block Diagram” window.

Next, move the cursor and **[Right-Click]** at the left node of the “Wait...” block. A menu should appear. Select **[Create]→[Constant]**, then change the value of the constant to: **10**. This means that the while loop will be activated after 10 milliseconds. (See Figure 19).

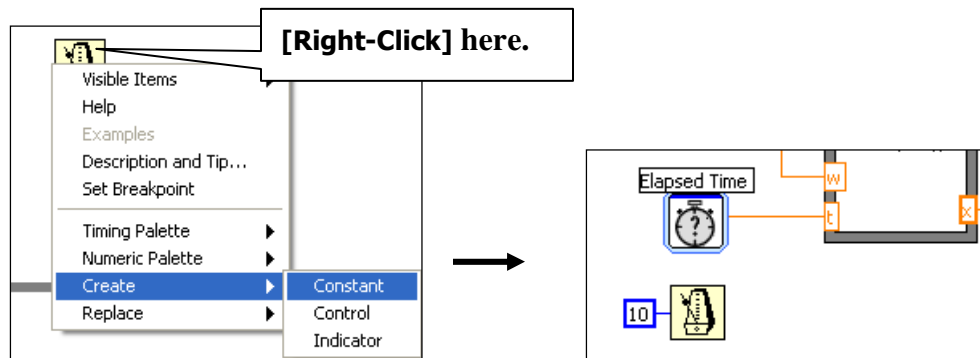


Figure 19. Attaching a constant value to the “Wait...” block.

8. Run/Test the VI project.

[Click] the “Run” button to run the VI (see Figure 20). Move the “Frequency” slider and observe what happens to the plot. You can **[Click]** the **[STOP]** button to terminate the run.

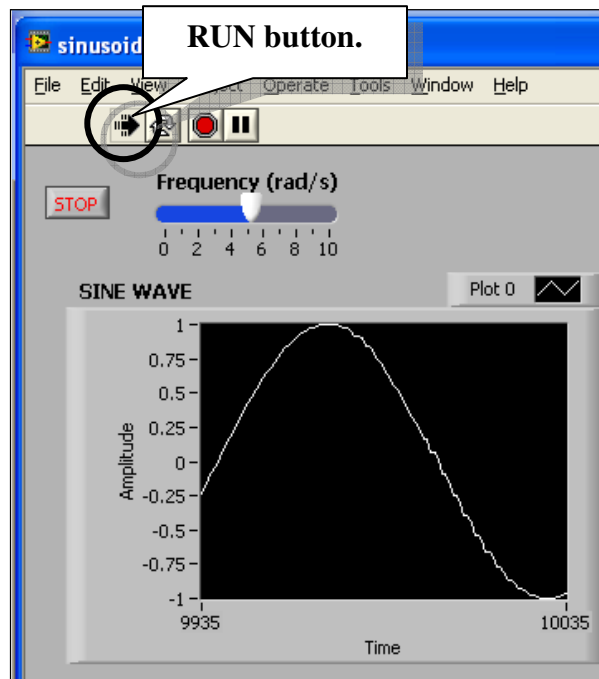


Figure 20. Running the VI project.

Remarks:

- i) The chart labels in the “SINE WAVE ” chart are the default labels. One usually needs to rename these labels.
- ii) The horizontal axis are in terms of “While loop” iterations. This means in our example, with the “Wait...” time of 10 milliseconds, the value of 9935 in Figure 20 means 99.35 seconds of elapsed time. Moreover, the range of samples shown in the chart is $(10035-9935)=100$ or 1 second.
- iii) It is advisable to save the VI frequently, specially when it works as desired.

9. Enhance/Modify the VI project and Add File Output.

- Include a numeric indicator to the “Frequency (slide)” control. (see Figure 21)
 - In the “Front Panel” window, access the Controls palette and expand to the **[Express]→[Numeric Indicators]** icon subdirectory. Select the **[Num Ind]** control and drop into the “Front Panel” window. Position the “Numeric Indicator” control to the right of the “Frequency (slide)” control.
 - **[Right-Click]** the “Numeric Indicator” control and select the **[Properties]** menu item, then match the entries given in Table 1.
 - In the “Block Diagram” window, make sure “Numeric Indicator” block is inside the “While loop” box. Position the cursor at a point in the wire connecting the “Frequency (slide)” block and the “Formula Node” block until the “wire spool” appears. Then drag a wire to connect it with the “Numeric Indicator” block.

Table 1. Properties used for
Numeric Indicator shown in Figure 21.

Tab	Item	Value
Display Format	Format	Floating Point
	Digits	3
	Precision Type	Digits of Precision
	Hide Trailing Zeros	<unchecked>
Appearance	Label : Visible	<unchecked>

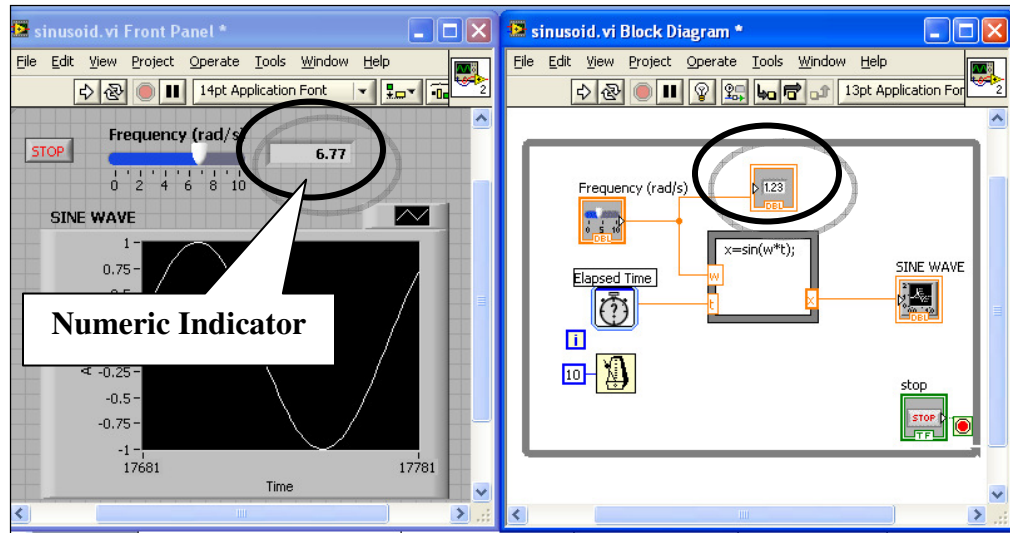


Figure 21. Adding a “Numeric Indicator” control to the “Frequency (slider)” control.

- Add a “Change to Dynamic Data” block. (see Figure 22)
 - In the “Block Diagram”, remove the wire connecting the “Function node” block and the “SINE WAVE” chart.
 - Access the “Functions palette” and expand the **[Express]→[Signal Manipulation]** icon subdirectory. Select **[to DDT]** icon (in some versions this is **[to Dynamic Data]**), then drag the block to somewhere between the “Function node” block and the “SINE WAVE” chart.
 - Wire the “**x**” output node to the input (left) node of the “to DDT” block. Also, wire the right (right) node of the “to DDT” block to the signal node of “SINE WAVE” chart. (Note that the second wire will be thicker. This type of wiring represents “dynamic data”, which means that time-information is now attached to the original data stream.)
 - **[Right-Click]** the “to DDT” block, then select the **[Properties]** menu item. Match the values given in Table 2.

Table 2. Properties used for
“to DDT” block in Figure 22.

<u>Item</u>	<u>Value</u>
Conversion	Single scalar
Start time	Now

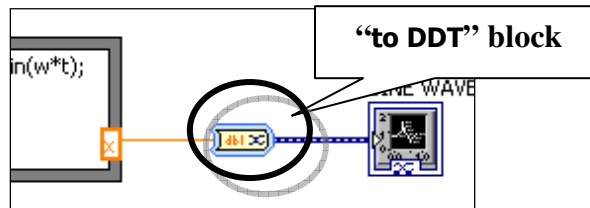


Figure 22. Insertion of a “to DDT” block which converts a single stream of data to dynamic data.

- Modify the properties of the “SINE WAVE” chart.
 - **[Right-Click]** the “SINE WAVE” chart in the “Front Panel” window, then select **[Properties]** menu item. Match the values given in Table 3. (see Figure 23)

Table 3. Modified Properties of “SINE WAVE” chart.

Tab	Variable	Item	Value
Format and Precision	Time(x-axis)	Type	Absolute time
		Time Format	Custom time format
		AM/PM	<on>
		HH:MM:SS	<on>
		Digits	1
		Date Format	Date unused
Scales	Time(x-axis)	Name	Time
		Autoscale	<unchecked>
		Minimum	0
		Maximum	10
	Amplitude (y-axis)	Name	x-values
		Autoscale	<unchecked>
		Minimum	-1
		Maximum	1

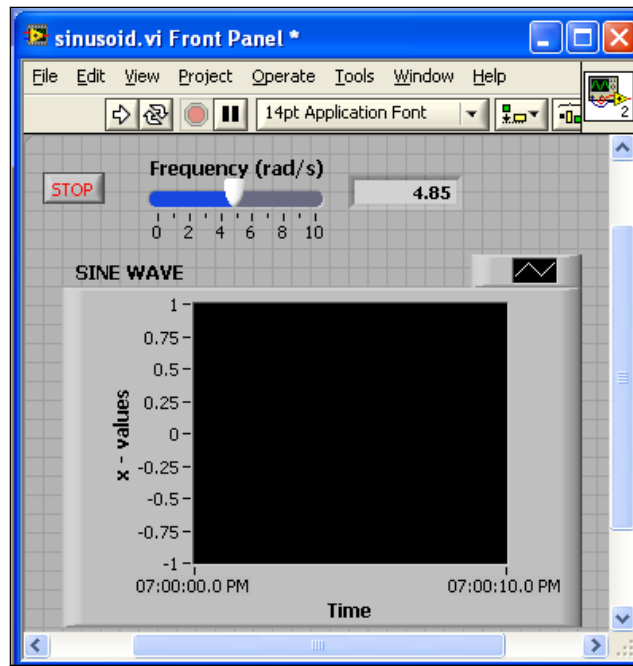


Figure 23. Modified chart properties.

- Add “Write Meas File” block (see Figure 24)
 - Access the “Functions palette” and expand the **[Express]→[Output]** icon subdirectory, then select the **[Write Meas File]** block and add inside the “While loop” box.
 - Reduce the block to an icon (**[Right-Click]** and select **[View as Icon]**) to save space.
 - Connect the “Signal” node of the “Write Meas File” block to the thin (orange) wire connecting “Function Node” block and the “to DDT” block.
 - **[Right-Click]** the “Write Meas File” block and select **[Properties]** menu item. Match the values given in Table 4.
 - Test the VI then check the output file using Excel. Note that the time included in the file will be the elapsed time, incremented by 0.01 seconds due to our chosen wait time of 10 milliseconds.

Table 4. Properties used for
“Write Meas File” block.

<u>Item</u>	<u>Value</u>
No headers	<on>
One column per channel	<on>

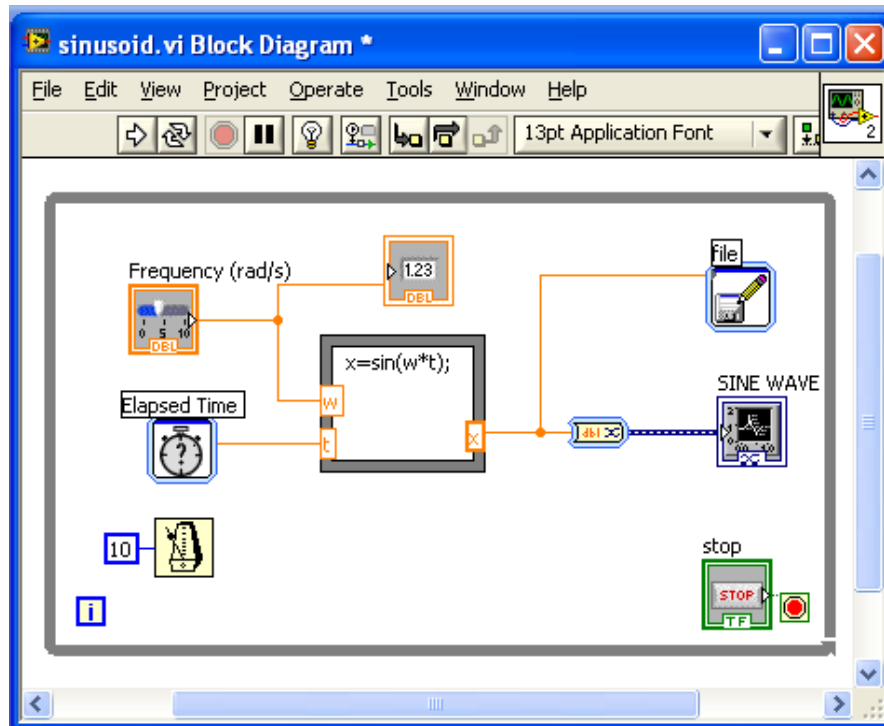


Figure 24. Final “Block-Diagram” window including a “Write Meas File” block.